



Viewing nCounter 360 Panel Web Reports


1. Save the .zip file to your desktop or a similar location on your computer.
2. Extract the contents by right-clicking on the .zip file and selecting **Extract All**, as shown below. Mac users may simply double click on the zipped file to initiate expansion.

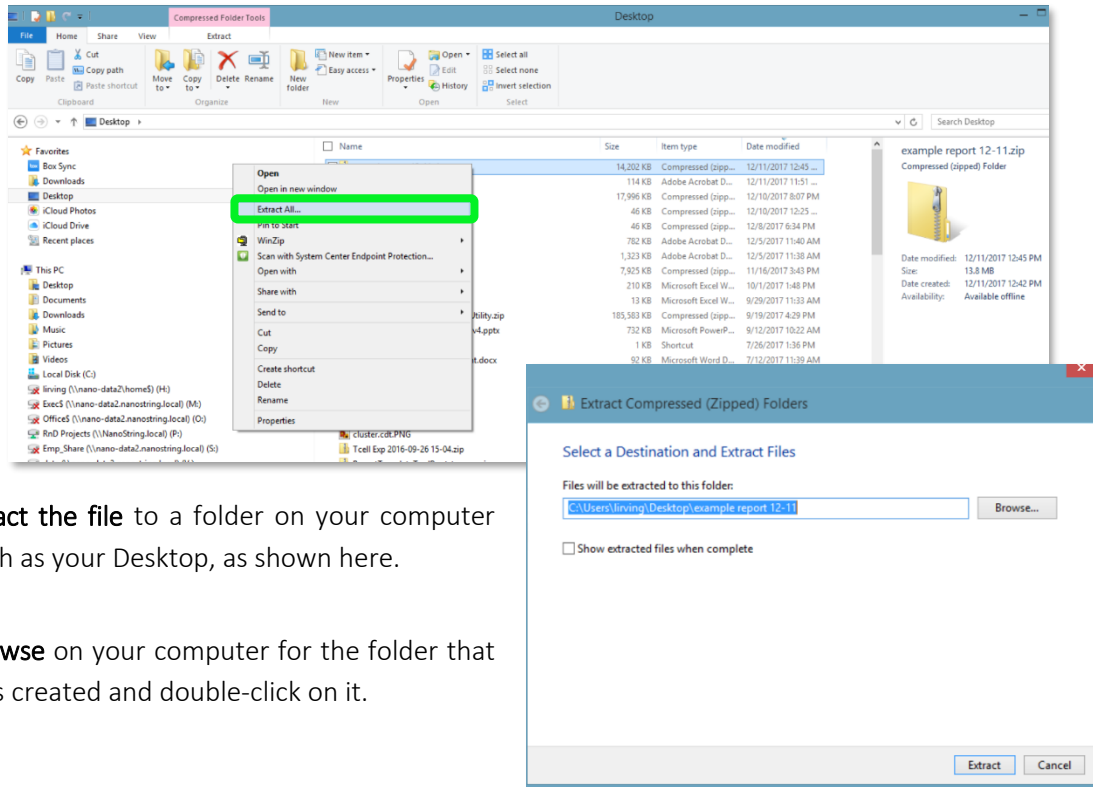


3. Extract the file to a folder on your computer such as your Desktop, as shown here.
4. Browse on your computer for the folder that was created and double-click on it.

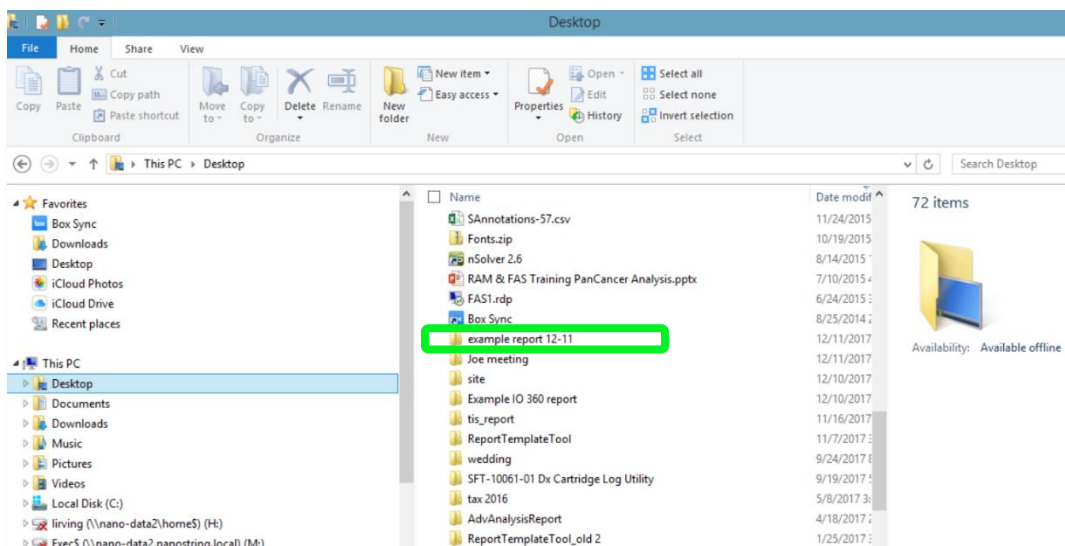


- Double-click on the **.html** document.
- Your browser should launch the report, as shown below. NOTE: Chrome is the recommended browser for this task.

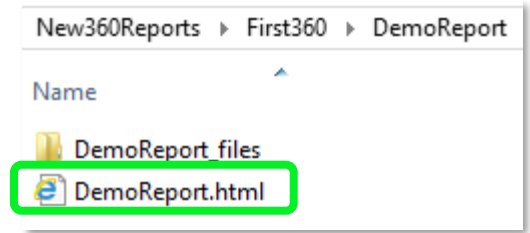

- Navigate** to different analyses using the panel menu on the left side of the window.

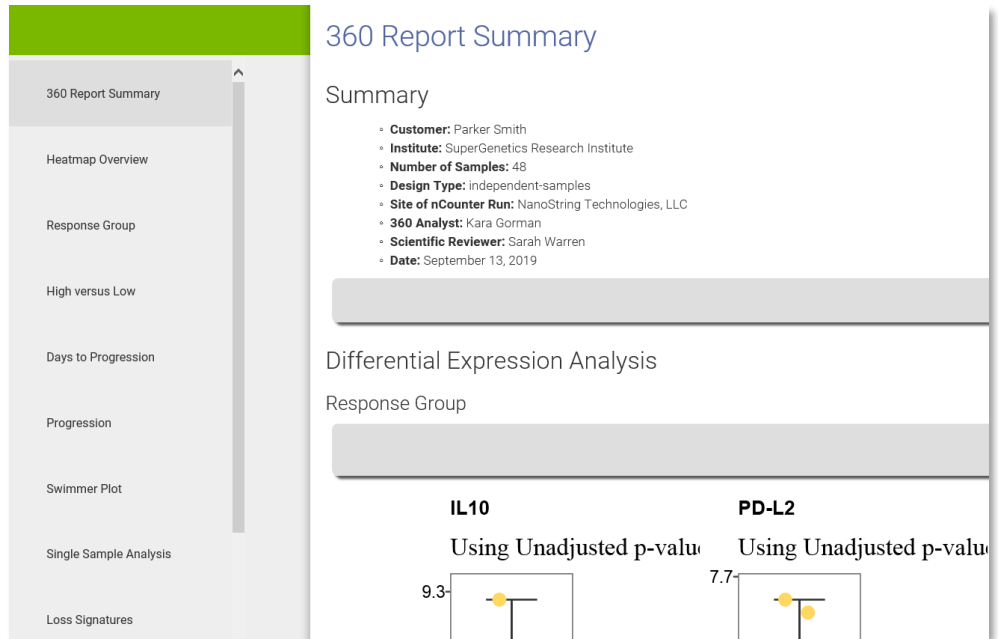

- Hover cursor** over upper right corner of any plot window to reveal a download button. Click to **download** a **.png** file of the present plot.

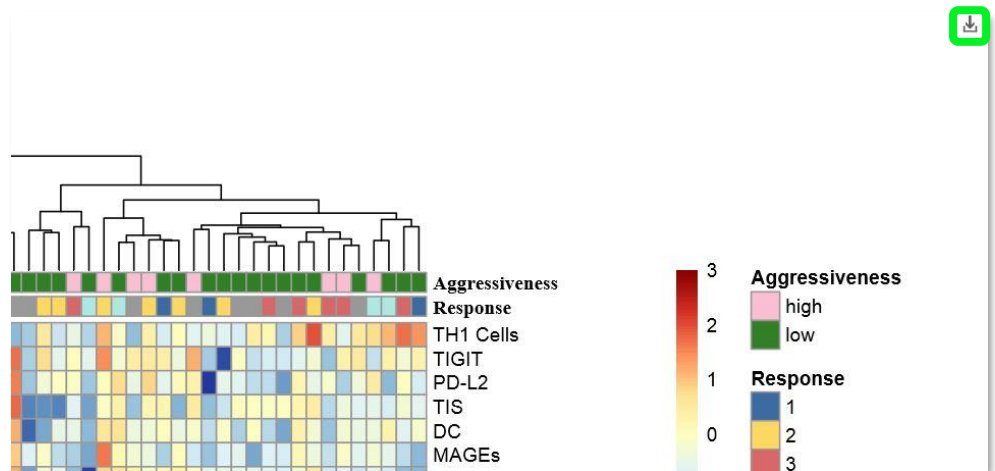

FOR RESEARCH USE ONLY. Not for use in diagnostic procedures

© 2019 NanoString Technologies, Inc. All rights reserved.

NanoString, NanoString Technologies, and the NanoString logo are trademarks or registered trademarks of NanoString Technologies, Inc., in the United States and/or other countries.